




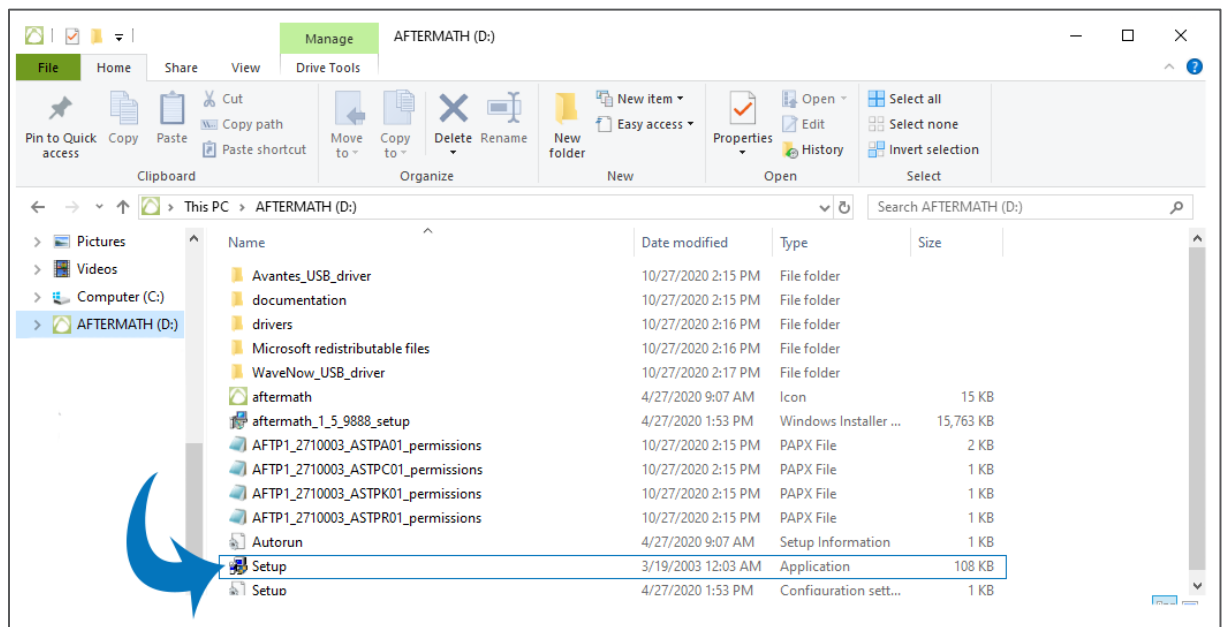



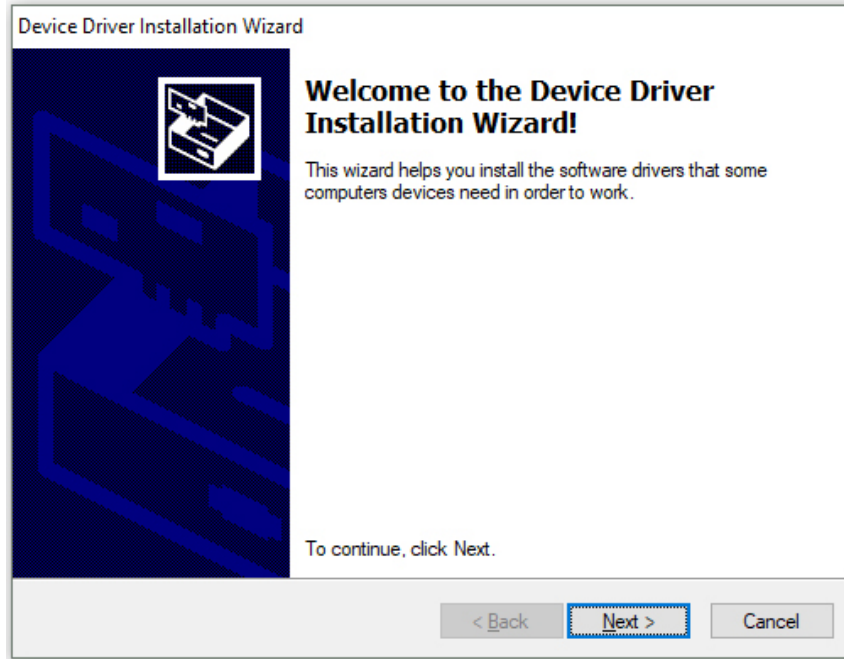


**READ ME.** Thank you for purchasing a great choice in potentiostats from Pine Research. We want your installation process to be fast and efficient. *It is essential to first install software and drivers before you connect the potentiostat to the computer.* Follow the Quick Start Installation instructions below.


-  **Find the User Guide.** This is typically a black binder located inside the main box or within the inner box if present.
-  **Leave the Potentiostat in the Box.** You will not need the instrument yet. Keep it safe until required in later steps.
-  **Obtain the USB Flash Drive.** It is located in a plastic flap inside the user guide. Connect it to your PC. 
-  **Install AfterMath Software.** Double-click the Windows Installer Package named "Setup" as shown below.






- 5  **Install Device Driver.** During the course of AfterMath installation, you will be prompted to install the device drivers.

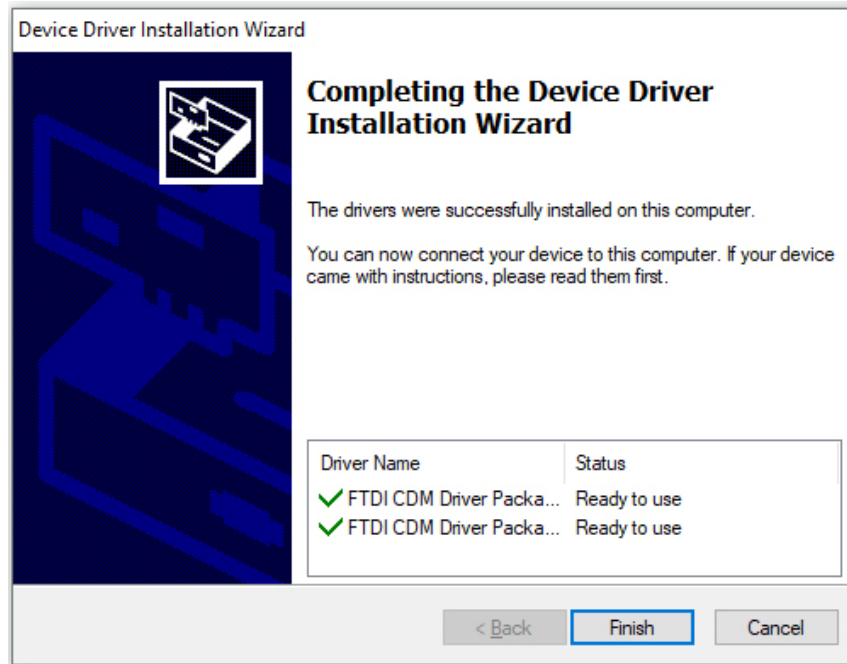


Verify the drivers were installed as shown:




Continue through any remaining steps and finish installation.

- 6  **Connect Potentiostat to the Computer by USB Cable.** Now that AfterMath software, instrument drivers, and support software have been installed, carefully unpack the potentiostat and connect it to your computer with the supplied USB cable.
  
- 7  **Power On Potentiostat.** Upon turning the potentiostat on, you should notice a Windows® notification (popup and/or sound) indicating that it detected new USB hardware. Wait about 30 seconds before proceeding to step 8.
  
- 8  **Start AfterMath.** You may have selected to place a shortcut on your desktop. If not, find where AfterMath was installed on your computer. Most commonly, the location is:  

**C:\Program Files (x86)\Pine\AfterMath**

 Open/start the aftermath.exe application in the folder.
  
- 9  **Verify the Status of the Potentiostat in AfterMath.** View the connected instruments in AfterMath. You should observe an entry for the potentiostat with its serial number in the *Idle* state.

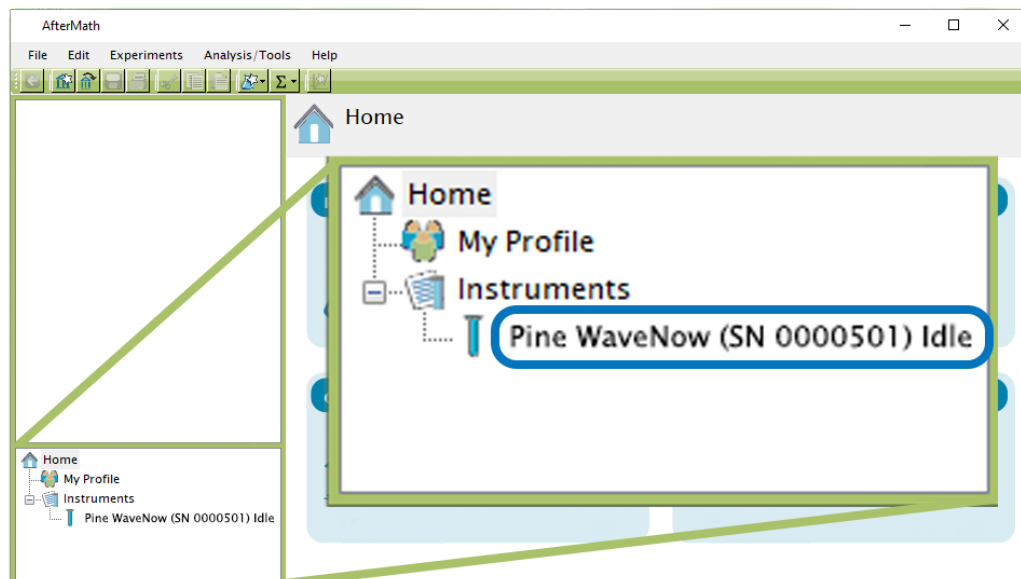

**Finished! If you have any issues or need assistance, please contact us**

 [pinewire@pineresearch.com](mailto:pinewire@pineresearch.com)

 **+1 (919) 782-8320**