



DO NOT BREAK EDGES WITH A FILE. USE ONLY 600 GRIT ABRASION SHEET BACKING BY A STRAIGHT EDGE.

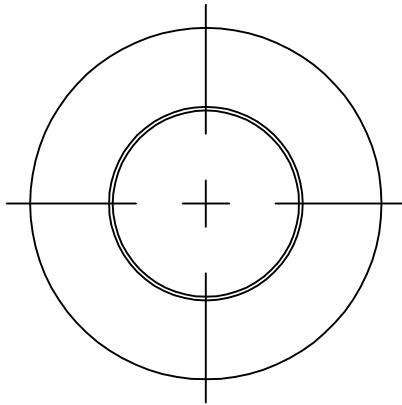

.005 x 45° CHAMFER BOTH SIDES

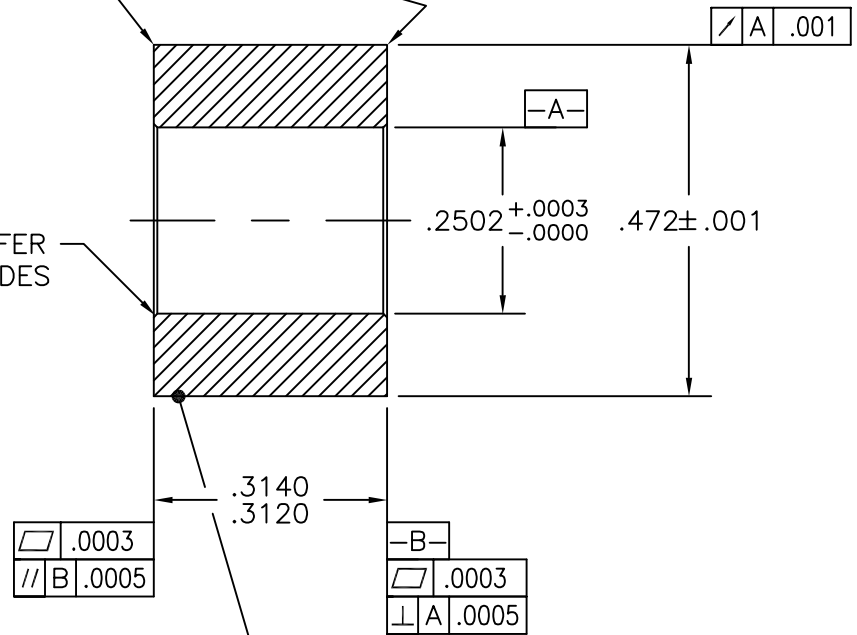

POLISH WITH 600 GRIT PAPER

NOTE: $\sqrt{16}$ ALL OVER.

TOLERANCES:	
X.XXX	= ±.005
X.XX	= ±.015
X.X	= ±.030
DEG.	= ±1/2°
FRAC.	= ±1/64
MACH. SURFACES $\sqrt{16}$ UNLESS NOTED	

SCALE: 4:1

2	5/17/93	REDRAWN	RSB
PINE		Instrument Company Grove City, PA. 16127-1091	
DWN.	RSB	CYLINDER 3 CM ² AREA	
CKD.	RAP		
ENG.	RAP	DWG. SIZE	DRAWING NO.
DATE	5/17/93	A	ACQC012CY
		SHEET	REV.
		1/1	2